




## Ductless Enclosures

- An Affordable, Effective Air Filtration System Solution
- Choose From Vertical or Low Profile Horizontal Design



Purair FUMEBOX-AP60V

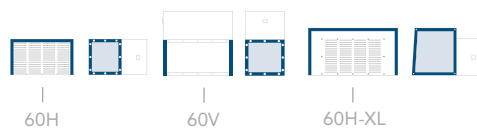

Schedule  
Contract # GS-47QSMS24D0042



## CONTENTS:

- Product Overview (p.2)
- Design Features (p.3)
- Performance & Selection (p.4)
- Containment & Filtration (p.5)
- Specifications (p.6)
- Options & Accessories (p.8)

## INTRODUCTION

Air Science® Purair® FUMEBOX ductless enclosures are designed to protect the user from chemicals, vapors or powders during low-volume chemical manipulations by effectively containing low concentrations of noxious fumes, vapors or powders.

## APPLICATIONS

Soldering \ Histology \ Cover Slipping \ Staining \ Pathology \ Graphic Arts, Sign Shops \ Art Conservation \ Electronics Assembly \ Particulate Control \ Manufacturing Processes \ Air Purification



*Deep into its second generation, Air Science embraces the diversity and cultural heritage of the founders and co-workers who are continuing a tradition of excellence. Demonstrating a commitment to adaptation, inclusion and quality output from a United States-based company with a domestic and global reach.*

## KEY FEATURES

- Low-profile, completely portable, no ductwork required.
- Compact size, easily positioned on shallow countertops or counters with overhanging wall cabinets.
- Clear viewing enclosure prevents chemical splash.
- Modular design permits multiple configurations.

## DUCTLESS TECHNOLOGY

### The Eco-Friendly Choice

Advanced carbon filtration technology offers a safe, high performance alternative to conventional ducted fume hoods for a broad range of applications.

- **Environmental Benefits.** Air Science ductless fume hoods isolate and trap chemical vapors to prevent ecological impact through release into the environment.
- **Versatile.** Each filtration system is selected for its specific application. Carbon filters are available in more than 14 configurations for use with vapors of organic solvents, acids, mercury and formaldehyde. HEPA/ULPA filters can be added.
- **Easy to Install.** The ductless fume hood is self-contained and does not require venting to the outside. Many units are portable and may be moved with minimal downtime and without filter changes. Set-up, operation and filter maintenance are straightforward.
- **Energy Efficient.** Because filtered air is returned to the room, no demands are required of the facility HVAC capacity for make-up air.
- **Cost Effective.** Facility ductwork, HVAC and construction costs are eliminated.
- **Safe to Use.** Cabinet airflow and face velocity protect users from incidental exposures to fumes.
- **Self-Testing.** (select models) Electronic airflow monitoring assures continuous safety. An electronic gas sensor monitors carbon filter performance.



Purair FUMEBOX-AP60H-XL

## CONTENTS:

- Product Overview (p.2)
- Design Features (p.3)
- Performance & Selection (p.4)
- Containment & Filtration (p.5)
- Specifications (p.6)
- Options & Accessories (p.8)



Purair FUMEBOX-AP60H

## DESIGN FEATURES

- A. Main On/Off Switch:** High quality rocker On/Off switch controls unit power.
- B. Steel Support Frame:** The chemical resistant epoxy coated steel frame adds mechanical strength.
- C. Electrostatic Pre-Filter:** The electrostatic pre-filter is accessible from inside the chamber and 91% effective down to 1-3 microns.
- D. Color:** The cabinet is white with blue trim; side and back panels are clear.
- E. Enclosed Filtration Chamber:** Main filters can be carbon and/or HEPA/ULPA to fit a variety of containment needs.

## ADDITIONAL FEATURES

**360 Degree Visibility:** Clear back and side panels allow ambient illumination into the chamber and provide users with an unobstructed view of its contents.

**Construction:** All models are available in polypropylene construction. See selection chart for specifications and dimensions. Available in 120V, 60Hz and 230V, 50Hz models.

## CONTENTS:

- Product Overview (p.2)
- Design Features (p.3)
- Performance & Selection (p.4)
- Containment & Filtration (p.5)
- Specifications (p.6)
- Options & Accessories (p.8)

Each Air Science fume hood includes features expressed through sound design and certified quality construction. Options and accessories add functional performance to meet specific applications.

## PERFORMANCE

The Air Science Multiplex™ Filter offers a range of options for high performance protection.


- Multiplex filter configuration permits a customized combination of filter media for a broad range of chemical families and biological agents if required.

A high capacity air handling system delivers face velocity of 100 fpm.

## DESIGN

Professional quality Air Science fume hoods comply with current technical and safety regulations. The cabinet frame and work surfaces, comprised of industrial components, are durable and chemically resistant.

The Air Science filter assembly is easy to access, easy to change, plus a unique filter clamping design eliminates bypass leakage outside the cabinet.

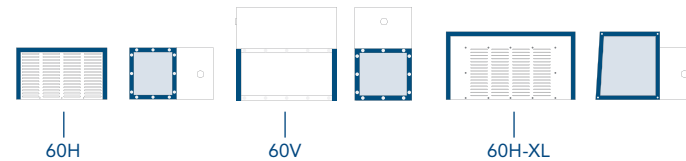

## SELECTION

Purair FUMEBOX ductless enclosures are available in 3 standard models, in metal or polypropylene construction, totaling 6 standard models.

## CONTROL

The **basic control panel** panel is standard and includes an On/Off switch and Filter Blockage alarm.



Basic Control Panel



Purair FUMEBOX-AP60V



The optional source capture hose converts the Purair FUMEBOX to a fume extraction unit with a 6" diameter (152 mm) hose to capture pollutants at their source. Hose can attach to any vented enclosure, exhaust port on your equipment or desk mount scoop.

## CONTENTS:

- Product Overview (p.2)
- Design Features (p.3)
- Performance & Selection (p.4)
- Containment & Filtration (p.5)
- Specifications (p.6)
- Options & Accessories (p.8)



## FILTRATION

The Purair FUMEBOX utilizes innovative filtration technology with the exclusive **Multiplex Filtration System**. The Multiplex Filtration System consists of a pre-filter, main activated carbon or HEPA/ULPA filter and safety activated carbon or HEPA/ULPA filter. The system permits a customized combination of filter media and configuration for chemical and physical adsorption specific to each application need.

The Air Science carbon filtration technique is based on enhanced, activated carbon particle formulations from specially selected, naturally occurring raw material that is superior to wood or other organic sources. The carbon is treated to attain the proper porosity and aggregate surface area and to react with several ranges of aerosolized chemicals moved through the filter by an air handling blower.

View available filters and descriptions on [page 7](#).

## FILTER CONFIGURATION

The Multiplex feature permits one or more filtration options to be combined to meet a wider range of multiple-use applications.

The Purair FUMEBOX can be equipped with a single activated carbon main filter. For safety against particulates, a HEPA or ULPA can be installed.



Filter disposal services are available in selected markets providing responsible destruction or recycling of saturated filters in authorized facilities.

The carbon filter is sized to fit the specified product model number and configured to optimize airflow across 100% of the filter surface area. The self-contained assembly maximizes filter efficiency, prolongs filter life, optimizes diffusion and saturation and improves user safety.

- P. Electrostatic Pre-Filter:** Protects the main filters from aerosols, mists, dust and particulates.
- C. Activated Carbon Main Filter:** A single filter configuration.
- H. HEPA/ULPA Filter, Optional:** Both HEPA and ULPA filters use micro-glass fiber media designed to capture fine particles and biologicals. Both filters can capture particles smaller than the micron size for which they are tested. HEPA and ULPA filter efficiencies are 99.97% at 0.3 microns and 99.999% at 0.12 microns respectively.

### MULTIPLY FILTRATION SYSTEM, SUMMARY

Application	Chemical	Powder/Biological
Primary Filter	C	H
Pre-Filter	P	P


The system can be configured for the capture of acids, bases and particulates, such as biological aerosols, when paired with HEPA or ULPA filters.

## AIRFLOW

The Purair FUMEBOX ductless enclosures maintain a constant face velocity of 100 fpm in compliance with USA and international standards for safety and performance. Contaminated air is pulled through the Multiplex filtration system; clean air is returned to the room.

The main filters are easy to replace and install. The filter clamps tightly against the filter gasket to prevent filter bypass and maintain filter integrity.


⚠ The pre-filter may be replaced while unit is in operation.

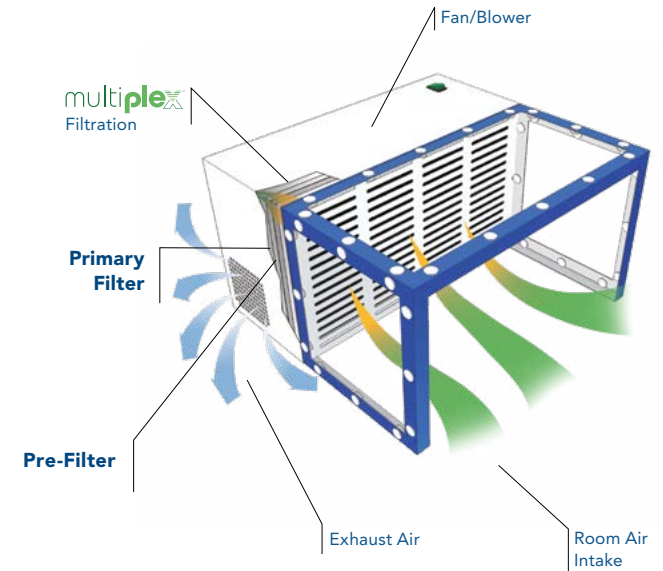

Through our partner company [Filtco Filters](#), Air Science is a single source supplier of all pre-filters, carbon filters and HEPA/ULPA filters used in our products.

**CONTENTS:**

- Product Overview (p.2)
- Design Features (p.3)
- Performance & Selection (p.4)
- Containment & Filtration (p.5)
- Specifications (p.6)
- Options & Accessories (p.8)

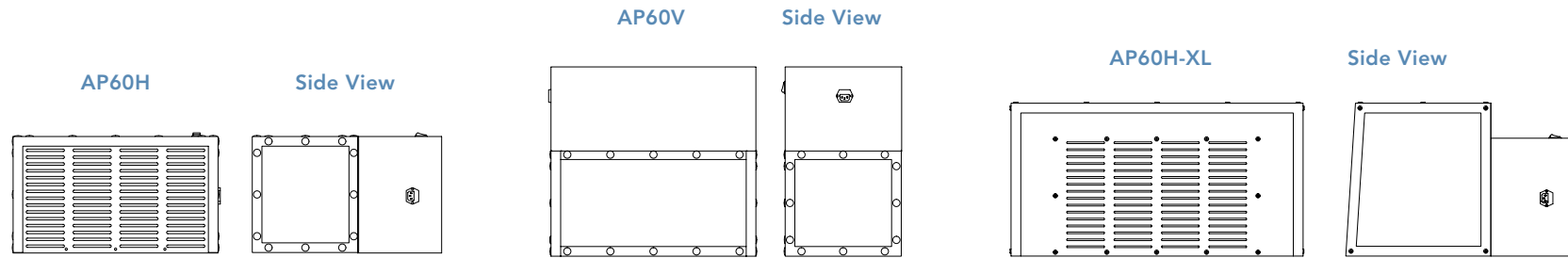

MODEL					WEIGHT (LBS/KG)	
Metal	Polypropylene	Internal (W x D x H)	External (W x D x H)	Shipping (W x D x H)	Net	Ship
AP60H	AP60H-PP	11.00" x 11.69" x 20.63" / 279 x 297 x 524 mm	20.72" x 12.02" x 21.16" / 526 x 305 x 537 mm	27" x 27" x 16" / 686 x 686 x 406 mm	45 / 20	60 / 27
AP60V	AP60V-PP	20.63" x 11.69" x 11.00" / 524 x 297 x 279 mm	21.16" x 12.02" x 20.72" / 537 x 305 x 526 mm	27" x 27" x 16" / 686 x 686 x 406 mm	45 / 20	60 / 27
AP60H-XL	AP60H-XL-PP	29.22" x 13.82" x 15.32" / 742 x 351 x 389 mm	29.75" x 24.72" x 15.67" / 756 x 628 x 398 mm	40" x 40" x 36" / 1016 x 1016 x 914 mm	50 / 23	100 / 45

**CONTENTS:**

- Product Overview (p.2)
- Design Features (p.3)
- Performance & Selection (p.4)
- Containment & Filtration (p.5)
- Specifications (p.6)
- Options & Accessories (p.8)

**PRODUCT SPECIFICATIONS**

<b>Filtration</b>	<b>AP60H</b>	<b>AP60V</b>	<b>AP60H-XL</b>
Airflow	<... 135.9 cfm ...>		
Face Velocity	<... 100 fpm ...>		
<b>Construction</b>	<b>AP60H</b>	<b>AP60V</b>	<b>AP60H-XL</b>
Finish	<... White epoxy-coated steel filtration unit with blue enclosure. Clear sides and top panels. ...>		
Blower	<... AC blower. ...>		
Controls	<... Main On/Off. ...>		
Electrical	<... 120V, 60Hz or 230V, 50Hz voltages available. Specify when ordering. Other voltage options available. ...>		
Monitoring	<... Filter blockage alarm, standard. ...>		
Pre-Filter Weight	<... 1 lb / 0.4 kg. ...>		
Filter Weight	<... 7 lbs / 3.5 kg. ...>		

**FILTER SPECIFICATIONS**

<b>Purair FUMEBOX</b>	<b>AP60H</b>	<b>AP60V</b>	<b>AP60H-XL</b>
Primary Filter(s)*	(1)	(1)	(1)
Pre-Filter*	(1)	(1)	(1)

\* For specific examples refer to Multiplex filtration system summary on [page 5](#).

**FILTER SUMMARY\***

<b>Formula</b>	<b>Description</b>
GP Plus!	The most widely used filter in the range, primarily for solvent, organic and alcohol removal.
ACI Plus!/ SUL	Designed to neutralize volatile inorganic acid vapors.
ACR	Iodine and methyl iodide vapors; It is frequently used for iodination reactions with lower level radioactive iodine.
ACM	Mercury vapor.
AMM	Removes vapors from dilute ammonia solutions and to remove low molecular weight amines.
FOR	Designed to oxidize formaldehyde and glutaraldehyde fumes; It is widely used in hospital pathology laboratories.
HEPA/UPLA	Powders and particulates.

\*Other formulas may be available.



## CONTENTS:

- Product Overview (p.2)
- Design Features (p.3)
- Performance & Selection (p.4)
- Containment & Filtration (p.5)
- Specifications (p.6)
- Options & Accessories (p.8)

## OPTIONS AND ACCESSORIES

Purair FUMEBOX		AP60H	AP60V	AP60H-XL
Spill Tray (Polypropylene)	Removable for easy cleaning.	TRAY-AP60	TRAY-AP60	TRAY-AP60-XL
Fume Extraction Unit		AP60EX		
Source Capture Hose*	Converts the Purair FUMEBOX to a fume extraction unit with a 6" diameter (152 mm) hose to capture pollutants at their source. Hose can attach to any, exhaust port on your equipment or desk mount scoop.		AP60EX	

\*Factory installed; specify when ordering.

## STANDARDS & COMPLIANCE

Quality Management Systems	ISO 9001:2015
Electrical Safety	UL-C-61010-1 CAN/CSA C22.2 61010-1-12 EN 61010-1:2010 CE Mark
OSHA, Occupational Safety and Health Information	OSHA Standard -29 CFR, Safety and Health Regulations for General Industry, 1910.1450: Occupational exposure to hazardous chemicals in laboratories. Part B, definition, laboratory type hood. This product may assist you with compliance or as part of your chemical hygiene plan. Please consult your Safety Officer and/or Industrial Hygienist.
Environment	ISO 14001:2015 ENERGY STAR® Partner



**CONTENTS:**

- Product Overview (p.2)
- Design Features (p.3)
- Performance & Selection (p.4)
- Containment & Filtration (p.5)
- Specifications (p.6)
- Options & Accessories (p.8)

**WARRANTY**

This product is protected by the Air Science Legacy Limited Lifetime Warranty™.



For details visit the [Warranty section](#) of our website.



120 6th Street \ Fort Myers, FL 33907  
T. 239-489-0024 \ Toll Free. 800-306-0656 \ F. 800-306-0677  
[www.airscience.com](http://www.airscience.com)

The information contained in this manual and the accompanying product are copyrighted and all rights are reserved by Air Science. Air Science reserves the right to make periodic minor design changes without obligation to notify any person or entity of such change.



Schedule  
Contract # GS-47QSMS24D0042

