




Compact Ductless Fume Adsorbers

- An Affordable, Effective Air Filtration System Solution
- Choose From Vertical or Low Profile Horizontal Design



Purair FUMEBOX-AP60V

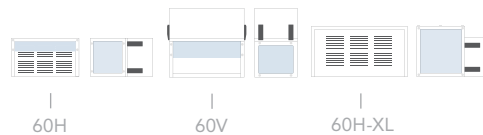

CONTENTS:

- Product Overview (p.2)
- Design Features (p.3)
- Performance & Selection (p.4)
- Containment & Filtration (p.5)
- Specifications (p.6)
- Options & Accessories (p.8)

Get a Quote. 

INTRODUCTION

Air Science® Purair® FUMEBOX compact ductless fume adsorbers are designed for nuisance vapor control, light-duty chemical handling and low-volume adsorption applications where portability, energy efficiency and filtration flexibility are desired.

APPLICATIONS

Soldering \ Graphic Arts, Sign Shops
Art Conservation \ Electronics Assembly
Particulate Control \ Manufacturing
Processes \ Air Purification



Deep into its second generation, Air Science embraces the diversity and cultural heritage of the founders and co-workers who are continuing a tradition of excellence. Demonstrating a commitment to adaptation, inclusion and quality output from a United States-based company with a domestic and global reach.

Air Science® USA LLC \ 120 6th Street, Fort Myers, FL 33907
T. 239-489-0024 \ airscience.com

Air Science® Technologies Ltd. \ United Kingdom
T. 0151 526 2457 \ airscience.com/UK

KEY FEATURES

- Low-profile, completely portable, no ductwork required.
- Compact size, easily positioned on shallow countertops or counters with overhanging wall cabinets.
- Clear viewing enclosure prevents chemical splash.
- Modular design permits multiple configurations.

DUCTLESS TECHNOLOGY

The Eco-Friendly Choice

Advanced carbon filtration technology offers a safe, high performance alternative to conventional ducted fume hoods for a broad range of applications.

- **Environmental Benefits.** Air Science ductless fume adsorbers isolate and trap chemical vapors to prevent ecological impact through release into the environment.
- **Versatile.** Each filtration system is selected for its specific application. Carbon filter cartridges are available in more than 14 configurations for use with vapors of organic solvents, acids and formaldehyde. HEPA/ULPA filter cartridges can be added.
- **Easy to Install.** The ductless fume adsorber is self-contained and does not require venting to the outside. Units are portable and may be moved with minimal downtime and without filter changes. Set-up, operation and filter maintenance are straightforward.
- **Energy Efficient.** Because filtered air is returned to the room, no demands are required of the facility HVAC capacity for make-up air.
- **Cost Effective.** Facility ductwork, HVAC and construction costs are eliminated.
- **Safe to Use.** Cabinet airflow and face velocity protect users from incidental exposures to fumes.



Purair FUMEBOX-AP60H-XL

Specifications are subject to change without notice or obligation on the part of Air Science. For questions contact Air Science.

CONTENTS:

- Product Overview (p.2)
- Design Features (p.3)
- Performance & Selection (p.4)
- Containment & Filtration (p.5)
- Specifications (p.6)
- Options & Accessories (p.8)

Get a Quote. 

PurairFUMEBOX™
Compact Ductless Fume Adsorbers

60H • 60V • 60H-XL
DESIGN FEATURES

3



Purair FUMEBOX-AP60H

DESIGN FEATURES

- A. Main On/Off Switch:** Simple to use on/off rocker switch controls unit power.
- B. Steel Support Frame:** The chemical resistant epoxy coated steel frame adds mechanical strength.
- C. Electrostatic Pre-Filter:** The electrostatic pre-filter cartridge is accessible from inside the chamber and 91% effective down to 1-3 microns.
- D. Color:** The cabinet is white; side and back panels are clear.
- E. Enclosed Filtration Chamber:** Main filter cartridge can be carbon or HEPA to fit a variety of containment needs.

ADDITIONAL FEATURES

High Visibility: Clear top and side panels allow ambient illumination into the chamber and provide users with an unobstructed view of its contents.

Construction: Available in 120V, 60Hz and 230V, 50Hz models.

CONTENTS:

- Product Overview (p.2)
- Design Features (p.3)
- Performance & Selection (p.4)
- Containment & Filtration (p.5)
- Specifications (p.6)
- Options & Accessories (p.8)

PurairFUMEBOX™
Compact Ductless Fume Adsorbers

60H • 60V • 60H-XL
PERFORMANCE & SELECTION

4

Get a Quote. 

Each Air Science ductless fume adsorber includes features expressed through sound design and certified quality construction. Options and accessories add functional performance to meet specific applications.

PERFORMANCE


The Air Science Multiplex™ Filtration System offers a range of options for high performance protection.

Multiplex filter configuration permits a customized combination of filter media for a broad range of chemical families and biological agents if required.

DESIGN

Professional quality Air Science ductless fume adsorbers comply with current technical and safety regulations. The cabinet frame and work surfaces, comprised of industrial components, are durable and chemically resistant.

The Air Science filter assembly is easy to access, easy to change, plus a unique filter clamping design eliminates bypass leakage outside the cabinet.

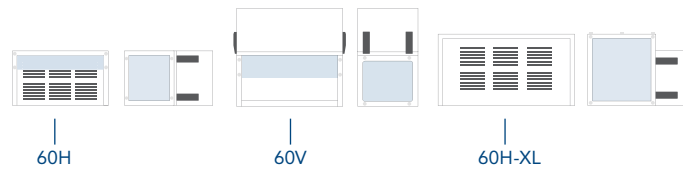

SELECTION

Purair FUMEBOX ductless fume adsorbers are available in 3 standard models, in metal construction, totaling 6 standard models.

CONTROL

The **basic control panel** is standard and includes an On/Off switch.



Purair FUMEBOX-AP60V



Basic Control Panel



The optional source capture hose converts the Purair FUMEBOX to a fume extraction unit with a 6" diameter (152 mm) hose to capture pollutants at their source. Hose can attach to any vented enclosure, exhaust port on your equipment or desk mount scoop.

CONTENTS:

- Product Overview (p.2)
- Design Features (p.3)
- Performance & Selection (p.4)
- Containment & Filtration (p.5)
- Specifications (p.6)
- Options & Accessories (p.8)

Get a Quote. 

multiple



FILTRATION

The Purair FUMEBOX utilizes innovative filtration technology with the exclusive **Multiplex Filtration System**. The Multiplex Filtration System consists of a pre-filter cartridge, main activated carbon or HEPA/ULPA filter cartridge and safety activated carbon or HEPA/ULPA filter cartridge. The system permits a customized combination of filter media and configuration for chemical and physical adsorption specific to each application need.

The Air Science carbon filtration technique is based on enhanced, activated carbon particle formulations from specially selected, naturally occurring raw material that is superior to wood or other organic sources. The carbon is treated to attain the proper porosity and aggregate surface area and to react with several ranges of aerosolized chemicals moved through the filter by an air handling blower.


View available filters and descriptions on [page 7](#).

FILTER CONFIGURATION

The Multiplex feature permits one or more filtration options to be combined to meet a wider range of multiple-use applications.

The Purair FUMEBOX can be equipped with a single activated carbon main filter cartridge. For safety against particulates, a HEPA or ULPA can be installed.

SECUR.
safe disposal service

 Filter disposal services are available in selected markets providing responsible destruction or recycling of saturated filters in authorized facilities.

The carbon filter cartridge is sized to fit the specified product model number and configured to optimize airflow across 100% of the filter surface area. The self-contained assembly maximizes filter efficiency, prolongs filter life, optimizes diffusion and saturation and improves user safety.

- P. Electrostatic Pre-Filter:** Protects the main filter cartridges from aerosols, mists, dust and particulates.
- C. Activated Carbon Main Filter:** A single filter cartridge configuration.
- H. HEPA/ULPA Filter, Optional:** Both HEPA and ULPA filter cartridges use micro-glass fiber media designed to capture fine particles and biologicals. Both filters can capture particles smaller than the micron size for which they are tested. HEPA and ULPA filter efficiencies are 99.97% at 0.3 microns and 99.999% at 0.12 microns respectively.

MULTIPLEX FILTRATION SYSTEM, SUMMARY


Application	Chemical	Powder/Biological
Primary Filter	C	H
Pre-Filter	P	P


The system can be configured for the capture of acids, bases and particulates, such as biological aerosols, when paired with HEPA or ULPA filters.

AIRFLOW

The Purair FUMEBOX ductless fume adsorber maintains a constant face velocity of 40 fpm in compliance with USA and international standards for safety and performance. Contaminated air is pulled through the Multiplex Filtration System; clean air is returned to the room.

The main filter cartridges are easy to replace and install. The filter clamps tightly against the filter gasket to prevent filter bypass and maintain filter integrity.

 The pre-filter cartridge may be replaced while unit is in operation.

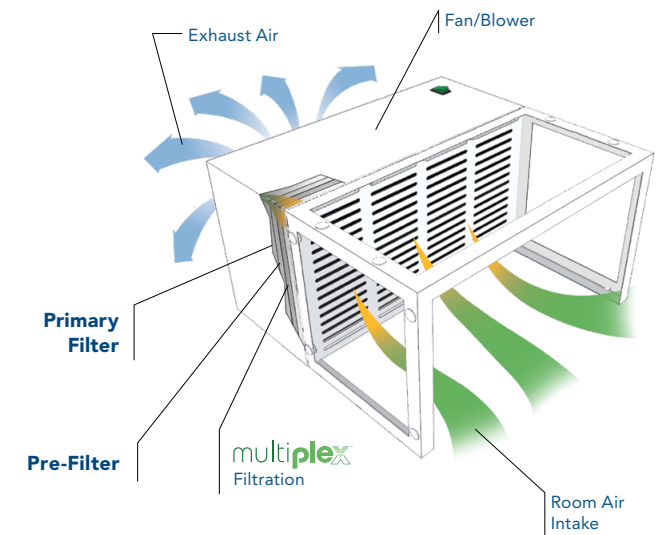


Through our partner company [Filtco Filters](#), Air Science is a single source supplier of all pre-filter cartridges, carbon filter cartridges and HEPA/ULPA filter cartridges used in our products.

Specifications are subject to change without notice or obligation on the part of Air Science. For questions contact Air Science.

CONTENTS:

- Product Overview (p.2)
- Design Features (p.3)
- Performance & Selection (p.4)
- Containment & Filtration (p.5)
- Specifications (p.6)
- Options & Accessories (p.8)

Get a Quote. 

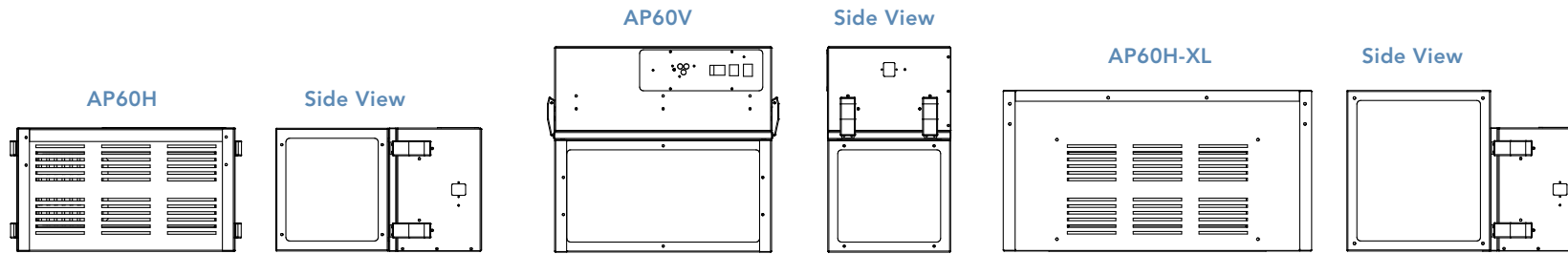

MODEL	VOLTAGE	DIMENSIONS			WEIGHT (LBS/KG)	
		Internal (W x D x H)	External (W x D x H)	Shipping (W x D x H)	Net	Ship
AP60H-A	120V, 60Hz	20.63" x 11.00" x 11.69" / 524 x 279 x 297 mm	20.72" x 12.02" x 21.16" / 526 x 305 x 537 mm	27" x 27" x 16" / 686 x 686 x 406 mm	45 / 20	60 / 27
AP60H-G	230V, 50Hz	20.63" x 11.00" x 11.69" / 524 x 279 x 297 mm	20.72" x 12.02" x 21.16" / 526 x 305 x 537 mm	27" x 27" x 16" / 686 x 686 x 406 mm	45 / 20	60 / 27
AP60V-A	120V, 60Hz	20.63" x 11.00" x 11.69" / 524 x 279 x 297 mm	21.16" x 12.02" x 20.72" / 537 x 305 x 526 mm	27" x 27" x 16" / 686 x 686 x 406 mm	45 / 20	60 / 27
AP60V-G	230V, 50Hz	20.63" x 11.00" x 11.69" / 524 x 279 x 297 mm	21.16" x 12.02" x 20.72" / 537 x 305 x 526 mm	27" x 27" x 16" / 686 x 686 x 406 mm	45 / 20	60 / 27
AP60H-XL-A	120V, 60Hz	30" x 13.82" x 15.32" / 742 x 351 x 389 mm	29.75" x 23.50" x 15.67" / 756 x 628 x 398 mm	40" x 40" x 36" / 1016 x 1016 x 914 mm	50 / 23	65 / 30
AP60H-XL-G	230V, 50Hz	30" x 13.82" x 15.32" / 742 x 351 x 389 mm	29.75" x 23.50" x 15.67" / 756 x 628 x 398 mm	40" x 40" x 36" / 1016 x 1016 x 914 mm	50 / 23	65 / 30

CONTENTS:

- Product Overview (p.2)
- Design Features (p.3)
- Performance & Selection (p.4)
- Containment & Filtration (p.5)
- Specifications (p.6)
- Options & Accessories (p.8)

Get a Quote. 

PRODUCT SPECIFICATIONS

Filtration	AP60H	AP60V	AP60H-XL
Airflow	41 cfm	41 cfm	62 cfm
Face Velocity	40 fpm / 0.2 m/s	40 fpm / 0.2 m/s	40 fpm / 0.2 m/s
Construction	AP60H	AP60V	AP60H-XL
Finish	<... White epoxy-coated steel filtration unit with blue enclosure. Clear sides and top panels. ...>		
Blower	<... EC blower. ...>		
Controls	<... Main On/Off. ...>		
Electrical	<... 120V, 60Hz or 230V, 50Hz voltages available. Specify when ordering. Other voltage options available. ...>		
Pre-Filter Weight	<... 1 lb / 0.4 kg. ...>		
Filter Weight	<... 7 lbs / 3.5 kg. ...>		
Sash Opening	<... 8" / 203 mm ...>		

FILTER SPECIFICATIONS

Purair FUMEBOX	AP60H	AP60V	AP60H-XL
Primary Filter(s)*	(1)	(1)	(1)
Pre-Filter*	(1)	(1)	(1)

* For specific examples refer to Multiplex filtration system summary on [page 5](#).

FILTER SUMMARY*

Formula	Description
GP Plus!	The most widely used filter in the range, primarily for solvent, organic and alcohol removal.
ACI Plus!/ SUL	Designed to neutralize volatile inorganic acid vapors.
AMM	Removes vapors from dilute ammonia solutions and to remove low molecular weight amines.
FOR	Designed to oxidize formaldehyde and glutaraldehyde fumes; It is widely used in hospital pathology laboratories.
HEPA/ULPA	Powders and particulates.

*Other formulas may be available.

CONTENTS:

- Product Overview (p.2)
- Design Features (p.3)
- Performance & Selection (p.4)
- Containment & Filtration (p.5)
- Specifications (p.6)
- Options & Accessories (p.8)

Get a Quote. 

OPTIONS AND ACCESSORIES

Purair FUMEBOX		AP60H	AP60V	AP60H-XL
Spill Tray (Polypropylene)	Removable for easy cleaning.	TRAY-AP60	TRAY-AP60	TRAY-AP60-XL
Fume Extraction Unit		AP60EX		
Source Capture Hose*	Converts the Purair FUMEBOX to a fume extraction unit with a 6" diameter (152 mm) hose to capture pollutants at their source. Hose can attach to any, exhaust port on your equipment or desk mount scoop.	AP60EX		

*Factory installed; specify when ordering.

STANDARDS & COMPLIANCE

Quality Management Systems	ISO 9001:2015
Electrical Safety	UL-C-61010-1 CAN/CSA C22.2 61010-1-12 EN 61010-1:2010 CE Mark
OSHA, Occupational Safety and Health Information	OSHA Standard -29 CFR, Safety and Health Regulations for General Industry, 1910.1450: Occupational exposure to hazardous chemicals in laboratories. Part B, definition, laboratory type hood. This product may assist you with compliance or as part of your chemical hygiene plan. Please consult your Safety Officer and/or Industrial Hygienist.
Environment	ISO 14001:2015 ENERGY STAR® Partner

CONTENTS:

- Product Overview (p.2)
- Design Features (p.3)
- Performance & Selection (p.4)
- Containment & Filtration (p.5)
- Specifications (p.6)
- Options & Accessories (p.8)

PurairFUMEBOX™
Compact Ductless Fume Adsorbers



60H • 60V • 60H-XL
OPTIONS & ACCESSORIES

Get a Quote.

Warranty Info.

WARRANTY

This product is protected by the Air Science Legacy Limited Lifetime Warranty™.



For details visit the [Warranty section](#) of our website.



120 6th Street \ Fort Myers, FL 33907
T. 239-489-0024 \ **Toll Free.** 800-306-0656 \ F. 800-306-0677
airscience.com

Air Science® Technologies Ltd. \ United Kingdom
T. 0151 526 2457 \ airscience.com/UK

The information contained in this manual and the accompanying product are copyrighted and all rights are reserved by Air Science. Air Science reserves the right to make periodic minor design changes without obligation to notify any person or entity of such change.



Schedule
Contract # GS-47QSMS24D0042

